

Data Sheet

Glucocorticoid Receptor Pathway GAL4 Reporter (Luc)-HEK293 cell Line **Catalog #: w70666**

Background

The glucocorticoid signaling pathway plays an important role in development, fluid homeostasis, cognition, immune response, and metabolism. Glucocorticoids are a class of steroid hormones that bind to the glucocorticoid receptor, causing it to translocate to the nucleus. Upon translocation, the receptor can regulate the transcription of a large number of genes, including those that regulate glucose metabolism and inflammatory responses.

Description

The *Glucocorticoid Receptor Pathway GAL4 Reporter (Luc) – HEK293 Cell Line* contains a firefly luciferase gene under the control of glucocorticoid receptor ligand binding domain that is fused to the DNA binding domain (DBD) of GAL4 (GAL4 DBD-GR) stably integrated into HEK293 cells. This fusion construct activates firefly luciferase expression under the control of a multimerized GAL4 upstream activation sequence (UAS). This allows for specific detection of glucocorticoid-induced activation of the glucocorticoid receptor without the need for individual transcriptional targets and with low cross-reactivity for other nuclear receptor pathways. This cell line is validated for response to stimulation of dexamethasone and to the treatment with mifepristone, an inhibitor of the glucocorticoid signaling pathway.

Applications

- Monitor glucocorticoid signaling pathway activity.
- Screen activators or inhibitors of the glucocorticoid signaling pathway.

Format

Each vial contains $\sim 1.5 \times 10^6$ cells in 1 ml of 10% DMSO.

Mycoplasma testing

The cell line has been screened using the PCR-based Venor™GeM Mycoplasma Detection kit (Sigma-Aldrich) to confirm the absence of *Mycoplasma* species.

Storage

Immediately upon receipt, store in liquid nitrogen.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.



9891 Irvine Center Dr. Suite 200
Irvine, CA 92618
United States
Tel: 1.800.831.1518
Fax: 1.800.831.1518
Email: sale@westbioscience.com

Culture conditions

Cells should be grown at 37°C with 5% CO₂ using MEM medium (Hyclone #SH30024.01) supplemented with 10% FBS (Life Technologies #26140-079), 1% non-essential amino acids (Hyclone #SH30238.01), 1 mM sodium pyruvate (Hyclone #SH30239.01), 1% Penicillin/Streptomycin (Hyclone #SV30010.01), 400 µg/ml of Geneticin (Life Technologies #11811031) and 50 µg/ml of Hygromycin B (Hyclone #SV3007001). If culturing cells in medium from other vendors, it may be required to lower the percentage of CO₂ in the incubator depending on the NaHCO₃ level in the basal medium.

It is recommended to quickly thaw the frozen cells from liquid nitrogen in a 37°C water-bath, and transfer to a tube containing 10 ml of growth medium without Geneticin or Hygromycin. Next, spin down the cells, resuspend cells in pre-warmed growth medium without Geneticin or Hygromycin, and transfer to a T25 flask. Culture the cells at 37°C in a 5% CO₂ incubator until the cells reach confluence. At first passage, switch to growth medium containing Geneticin and Hygromycin. Cells should be split before they reach complete confluence.

To passage the cells, rinse with phosphate-buffered saline (PBS) and detach cells from the culture vessel with Trypsin/EDTA. After detaching, add complete growth medium and transfer to a 15 ml conical tube to spin down the cells. After spinning, resuspend cells in complete growth medium and seed appropriate aliquots of cell suspension into new culture vessels.

Note: Just after thawing and at low density, the cells may grow at a slower rate. It is recommended to split the cells with ~1:4 ratio at the beginning of culturing. After several passages, the cell growth rate increases and the cells can be split with 1:8 -1:20 ratio weekly.

Functional Validation and Assay Performance

The following assays are designed for a 96-well format. To perform the assay in different tissue culture formats, the cell number and reagent volume should be scaled appropriately.

Materials Required but Not Supplied


- Dexamethasone (Sigma #D4902)
- Mifepristone/RU-486 (Sigma #M8046)
- HEK293 assay medium: DMEM (without phenol red) (Hyclone #SH30604.01) + 10% charcoal/dextran-treated FBS (Hyclone #SH30068.02), 4 mM L-Glutamine + 1% Pen/Strep
- 96-well tissue culture treated white clear-bottom assay plate (Corning #3610)
- One-Step Luciferase Assay System (West Bioscience Cat. #w70701)
- Luminometer

A. Dose response of *GAL4 Reporter (Luc)* – HEK293 cells to dexamethasone

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

1. Harvest GAL4 reporter (Luc)-HEK293 cells from culture in growth medium and seed cells at a density of ~30,000 cells per well into a white clear-bottom 96-well microplate in 45 μ l of assay medium.
2. Incubate cells at 37°C in a CO₂ incubator overnight (~16 hours).
3. Add 5 μ l of three-fold serial dilutions of dexamethasone in assay medium to stimulated wells. Typically, 8 – 12 concentrations are used.
Add 5 μ l of assay medium to the unstimulated control wells.
Add 50 μ l of assay medium to cell-free control wells (for determining background luminescence).
Set up each treatment in at least triplicate.
4. Incubate the plate at 37°C in a CO₂ incubator for 24 hours.
5. Perform luciferase assay using One-Step Luciferase Assay System and following the protocol provided: Add 100 μ l of One-Step Luciferase reagent per well and rock at room temperature for ~10 minutes. Measure luminescence using a luminometer. *If using other luciferase reagents from other vendors follow the manufacturer's assay protocol.*
6. Data Analysis: Subtract the average background luminescence (cell-free control wells) from the luminescence reading of all wells. Fold induction of GAL4 luciferase reporter expression = background-subtracted luminescence of dexamethasone-stimulated well / average background-subtracted luminescence of unstimulated control wells

Figure 1. Dose response of GAL4 reporter (luc)-HEK293 cells to dexamethasone. The results are shown as fold induction of GAL4 luciferase reporter expression.

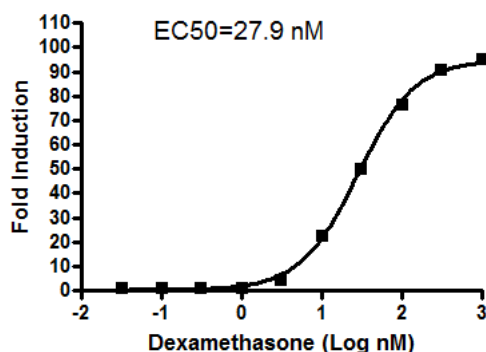


OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

B. Inhibition of dexamethasone-induced reporter activity by an antagonist of the glucocorticoid signaling pathway in *GAL4 Reporter (Luc)-HEK293 cells*

1. Harvest GAL4 reporter (Luc)-HEK293 cells from culture in Growth medium and seed cells at a density of ~30,000 cells per well into a white clear-bottom 96-well microplate in 45 μ l of assay medium.
2. Incubate cells at 37°C in a CO₂ incubator overnight (~16 hours).
3. Add 5 μ l of diluted mifepristone (a glucocorticoid antagonist) in three-fold serial dilutions in assay medium to inhibited wells. Typically 8 – 12 concentrations are used. Set up each treatment in at least triplicate. Incubate cells at 37°C in a CO₂ incubator for 1 hour.
4. Add 5 μ l of 200 nM dexamethasone in assay medium to stimulated wells (final dexamethasone concentration = 20 nM).
Add 5 μ l of assay medium to the unstimulated control wells (for determining the basal activity).
Add 55 μ l of assay medium to cell-free control wells (for determining background luminescence).
Set up each treatment in at least triplicate. Incubate the plate at 37°C in a CO₂ incubator for 24 hours.
5. Perform luciferase assay using ONE-Step Luciferase Assay System and following the protocol provided: Add 100 μ l of One-Step Luciferase reagent per well and rock at room temperature for ~10 minutes. Measure luminescence using a luminometer. *If using other luciferase reagents from other vendors follow the manufacturer's assay protocol.*
6. Data Analysis: Obtain the background-subtracted luminescence by subtracting the average background luminescence (cell-free control wells) from the luminescence reading of all wells.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

Figure 2. Inhibition dose response curve of dexamethasone-induced reporter activity by mifepristone in GAL4 reporter (Luc)-HEK293 cells. The results are shown as percentage of luminescence. The background-subtracted luminescence of cells stimulated with dexamethasone in the absence of mifepristone was set at 100%.

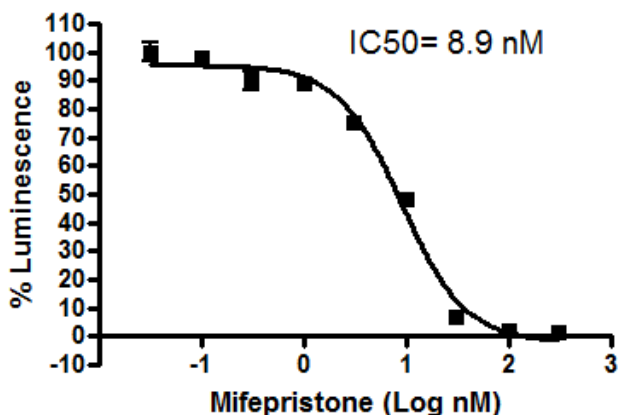

Reference

Paguio A, *et al.* (2010) Improved Dual-Luciferase Reporter Assays for Nuclear Receptors. *Curr Chem Genomics*. **4**: 43-49.

License Disclosure

Purchase of this cell line grants you with a 10-year license to use this cell line in your immediate laboratory, for research use only. This license does not permit you to share, distribute, sell, sublicense, or otherwise make the cell line available for use to other laboratories, departments, research institutions, hospitals, universities, or biotech companies. The license does not permit the use of this cell line in humans or for therapeutic or drug use. Inappropriate use or distribution of this cell line will result in revocation of the license and result in an immediate cease of sales and distribution of West products to your laboratory. Publications using this cell line should reference West Bioscience, Inc., San Diego.

OUR PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.